





Tutorial Guide 1 – Setup and Application Launch



This page is left intentionally blank

1. INSTALLATION

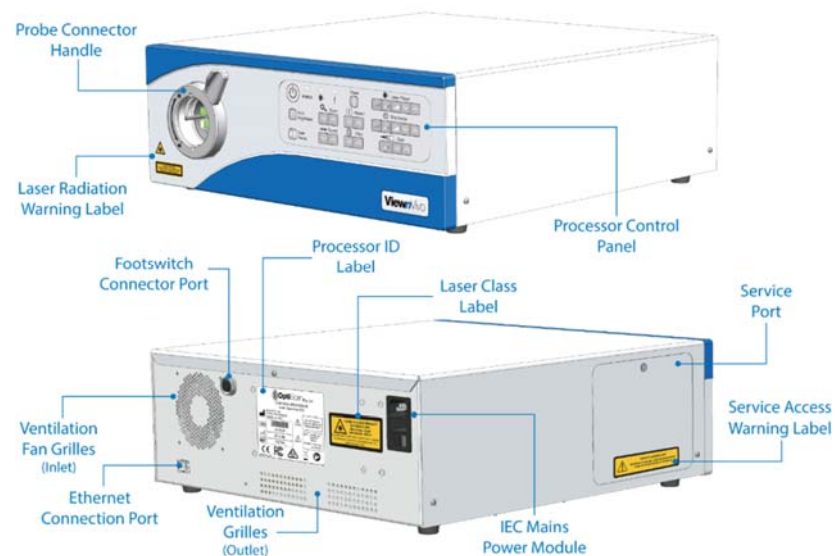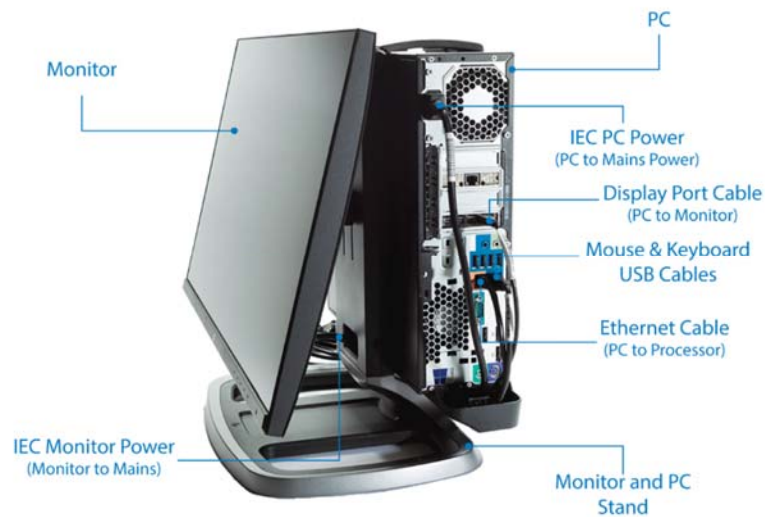

Connect PC Cables



Connect Viewvivo Processor Cables


2. POWER UP THE SYSTEM & CONNECTING THE PROBE



Turn the Viewvivo Processor Main Power ON




Turn on the PC

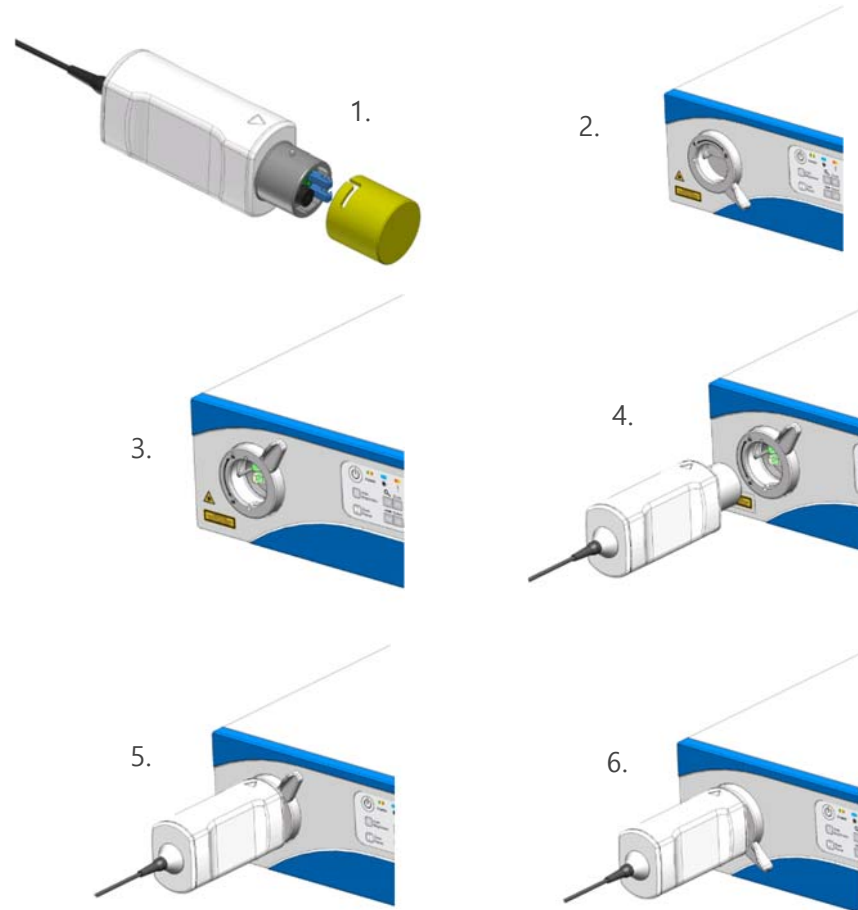


Connecting the Probe

Mains Power is OFF	Mains Power ON System Turned OFF	System Turned ON
<p>POWER</p>	<p>POWER</p>	<p>POWER</p>

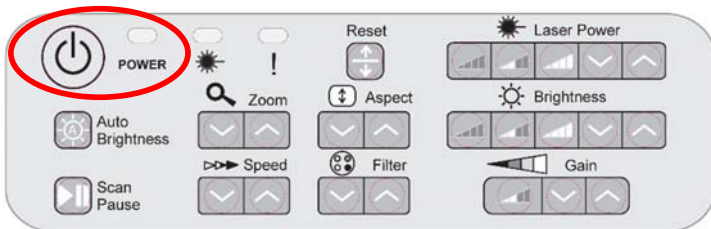

The Processor Control Panel

3. LAUNCHING THE IMAGER SOFTWARE

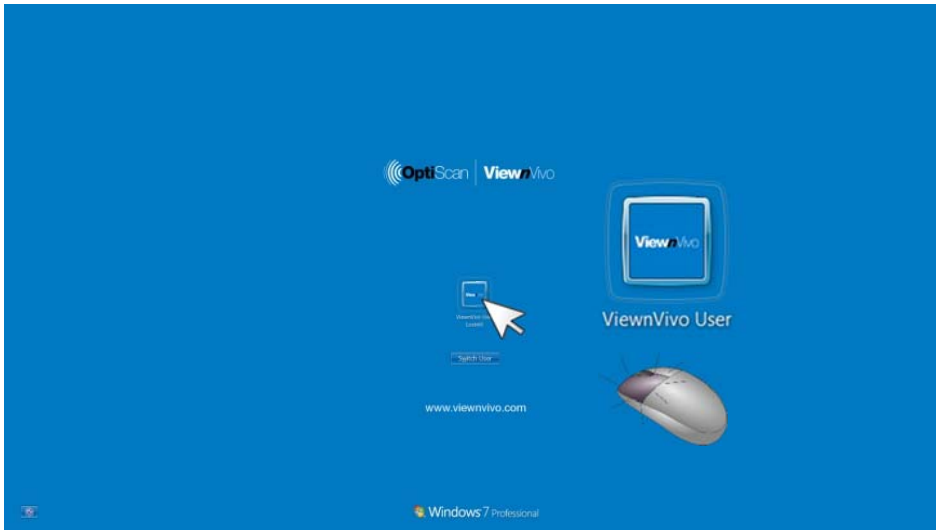

Step 1: Login using the 'ViewnVivo User' account – No Password is needed



Step 2: At the Windows OS Desktop, Double click on the Imer Application Icon



Step 3: The Application is trying to connect to the Confocal Processor



Step 4: The Application has established communications with the device

2. DEVICE WARMUP




Step 5: The System is Initialising

Note: If the 'Confocal Probe' is not connected a "No Endoscope" status message will appear in the bottom left hand corner, and a "LaserInterLockOpen" status message will appear on the top left-hand corner and the 'System Status LED' on the 'Processor Control Panel' will be coloured "Red". Connect a Probe to continue initialisation




No Endoscope

Status: LaserInterLockOpen 0 μ w

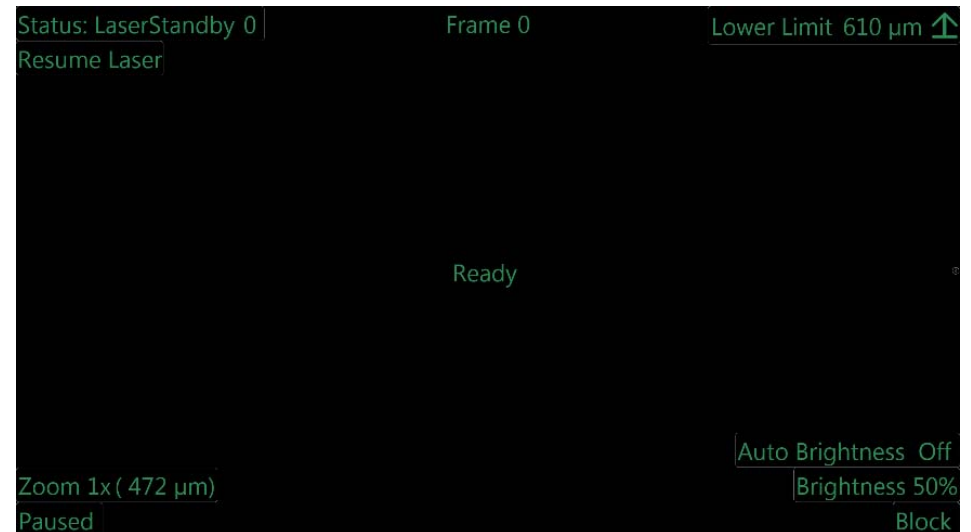

Step 6: The System is Ready to Use, and all preferences have been loaded

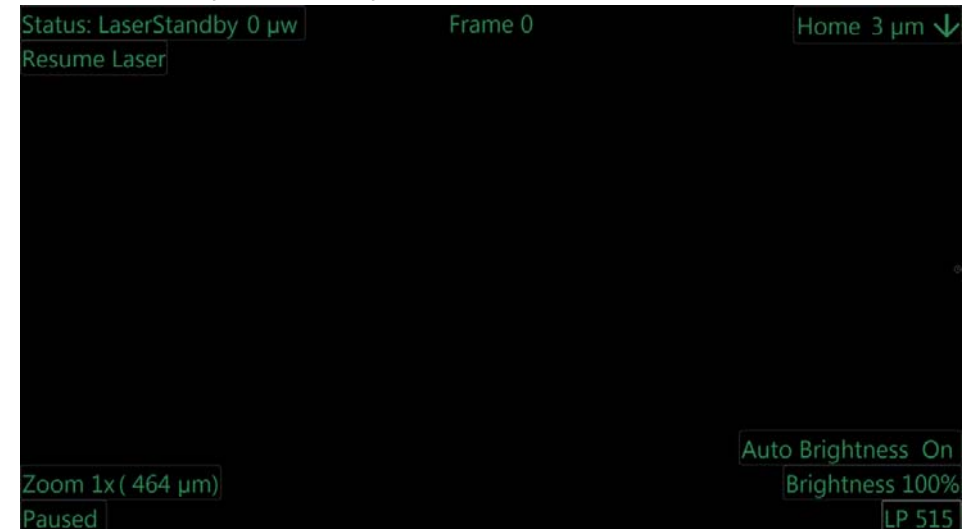

Step 7: The System is Ready to Use, and all preferences have been loaded

This page is left intentionally blank

Manufactured by



Optiscan Pty. Ltd.

Head Office

16 Miles Street, Mulgrave

Victoria 3170, AUSTRALIA

Phone: +61 3 9538 3333

Fax: +61 3 9562 7742

www.optiscan.com

Sales

sales@optiscan.com

Customer Support

service@optiscan.com

Additional Resources and Links

www.optiscan.com

www.viewnvivo.com